

RAZOR® EX and RAZOR® Mk II Battery Compatibility

Summary

This Technical Note is to advise RAZOR users about battery compatibility between instrument versions.

The RAZOR Mk II battery is compatible with the RAZOR EX, but it is not compatible with the RAZOR EX battery charger. Use the table below to identify which battery accessories are compatible.

RAZOR EX customers who need a new battery must order the new battery pack and a compatible charger.

Part Number	Description
RAZR-ASY-5100	RAZOR Mk II Battery Pack
RAZR-ASY-5200	RAZOR Mk II Battery Charger

RAZOR Battery Compatibility

The table below indicates which RAZOR battery accessories can be used with the RAZOR Mk II or the RAZOR EX.

	RAZOR Mk II Instrument	RAZOR Mk II Battery	RAZOR Mk II Charger
RAZOR EX Instrument	-	Compatible*	-
RAZOR EX Battery	Compatible	-	Not Compatible
RAZOR EX Charger	-	Not Compatible	-

***NOTE:** If using a RAZOR Mk II battery with a RAZOR EX instrument the user may experience an anomalous “voltage too high” warning. This is the result of newer and more efficient battery technology. This can be resolved by rebooting the instrument until the battery level drops to RAZOR EX starting levels.

TECHNICAL NOTE

Technical Support Contact Information

BioFire Defense is dedicated to providing the best customer support available. If you have any questions or concerns about this information, please contact our Technical Support team for assistance.

General Information

Email: support@biofiredefense.com

Phone: 1-801-262-3592



TECHNICAL
::: NOTE